


## ASSEMBLY INSTRUCTIONS FOR BUILT-IN FUME HOOD HI2P


1. Draw a rectangle with the dimensions of the HI2P hood (350mm X 425mm) on the table.
2. Make 4 holes with a hole saw or drill bit, then cut out the rectangle with a jigsaw.

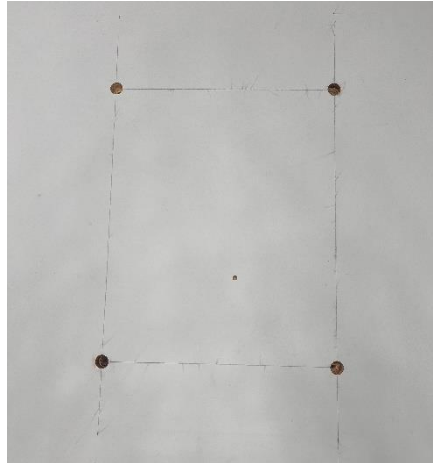

3. Insert the hood fitted with its metal corners, until it rests on the table. If necessary, adjust the hole dimensions with light grinding.
4. Locate the 4 grid fixing holes



5. Drill the 4 holes crossing the work surface with a diameter of 6mm.

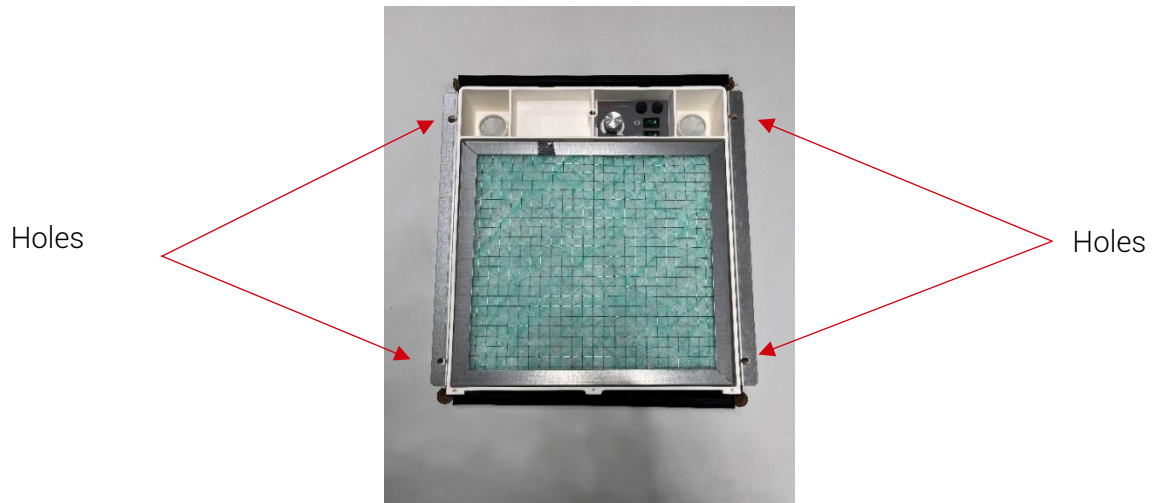

6. Install the filter and pre-filter (optional), then position the grid on the table until it stops. If necessary, complete the installation with 4 nuts to screw onto the grid studs (for table thickness 20 to 30mm).

