

## OPTION : MOTION DETECTION TRIGGERING

The **HI9P**, **HI5VP**, **HI7VP** exhaust hoods offer an optional automatic lighting and motor activation function.



HI5VP : straight fume hood



HI9P : overhead hood




HI7VP : pyramidal fume hood

### FOCUS ON THE FUME HOOD

Detector control button

Detection slot

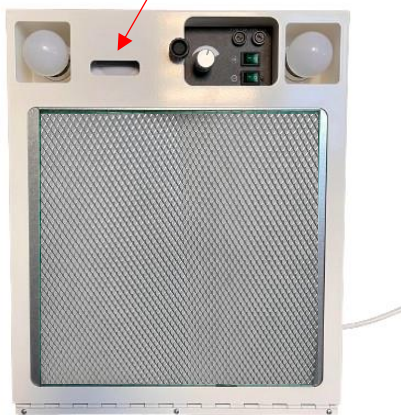

view with cache

Detector control button

Detection device



view without cache

## Working Principle:

- When movement is detected in front of the extraction box, the motor and the hood lighting turn on.
- The engine speed is established at the speed set by the potentiometer
- If movement is detected again during the active period, the activation delay is extended by a full period.
- We recommend setting the operating time to the maximum duration (> 5 minutes). To do this, turn the button on the left completely to the right. This setting guarantees good evacuation of particles and chemical agents for a few minutes after the end of handling.
- For information, the light sensitivity is adjusted beforehand in our workshops, so **no further manipulation is required.**

